

## MAGE3/MHC mRNA-LNP

Ready-to-use lipid nanoparticles

### Order Information

Catalog#	Size	GenPept No.
PM-LNP-0015	200uL	

### Description

Melanoma-associated antigen 3 (MAGE3 or MAGEA3), a protein belonging to the melanoma-associated antigen gene family, has been found in many tumors, including melanoma, non-small cell lung cancer, hematological malignancies, and others. MAGE genes are activated in tumors of many different histological types. Due to this strict tumor specificity, MAGE antigens are of particular interest in cancer immunotherapy. The major histocompatibility complex (MHC) is a large locus on vertebrate DNA that contains a tightly linked set of polymorphic genes that encode cell surface proteins critical to the adaptive immune system. Classical MHC molecules present epitopes to the TCR of CD8+ T lymphocytes. The antigens/ligands of many of these molecules are still unknown, but they can interact with

### Composition

mRNA-LNPs suspended in PBS (-Ca, -Mg) (pH: 7.0-7.4).

### Translated Protein sequence


Available up request

#### Products and Services

- Monoclonal Antibody
- Bispecific Antibody
- mRNA-LNP
- Hybridoma Sequencing
- CAR-T/NK
- Stable Cell Line

## **MAGE3/MHC mRNA-LNP**

### **Ready-to-use lipid nanoparticles**

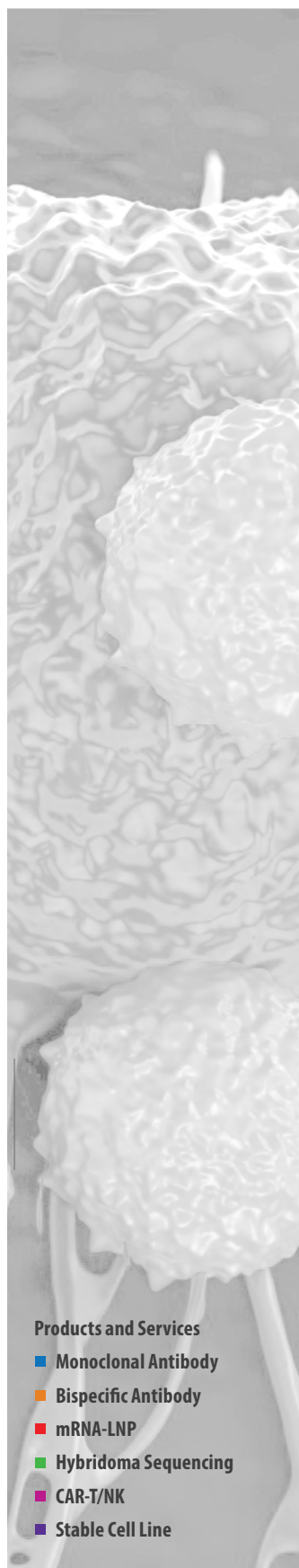

#### **Products and Services**

- **Monoclonal Antibody**
- **Bispecific Antibody**
- **mRNA-LNP**
- **Hybridoma Sequencing**
- **CAR-T/NK**
- **Stable Cell Line**

#### **Storage**

Product is delivered on blue ice. Store at 4°C for up to 3 months.

#### **Application & Handling**

Upon receiving product, briefly pulse spin before opening to ensure product is at bottom of container. It is important not to spin for too long as this may rupture mRNA-LNPs. Do not vortex. Work with mRNA-LNPs on ice and minimize the time that the product spends at room temperature. After handling the product during experiments, return immediately to ice. mRNA-LNP products should only be handled with certified RNase-free reagents and consumables. Use of filtered pipette tips is highly recommended.

#### **Safety & Research Disclosure**

All ProMab mRNA lipid nanoparticle products are for in vitro research use only. Products are not FDA approved for human use.

#### **General Protocol**

1. Prior to transfection: Plate 1ml of cells at a density of [1.0E6 cells/ml] in a single well of a 12-well culture plate. Ensure the cells you are using are viable and healthy. Try not to let your cells sit for longer than 5 minutes prior to transfection. Cell clumping at the time of transfection may reduce transfection efficiency. \*Note: If cell clumping occurs, gently pipette your culture up & down to ensure you have a single cell suspension before transfecting.
2. Briefly pipette mRNA-LNP mix up & down to resuspend. Add 20-40ul of the mRNA-LNP product dropwise directly to your 1ml culture. Gently tilt plate back and forth to mix (not necessary if you are using cells which will be immediately placed back into a shaker). Place your transfected cells back into their original culture conditions.
3. Check cell expression by FACS or by using other detection methods at 24hr intervals after transfection. \*Note: This is a generalized protocol for transfection using mammalian suspension culture cells. Transfection volume may be scaled up or down proportionately using the volumes given. HEK-293s cells were grown and transfected in FreeStyle™ F17 Expression Medium (Gibco, Cat#: A1383501), supplemented with GlutaMAX™ (Gibco, Cat#: 35050061), and Poloxamer 188 Non-ionic Surfactant (Gibco, Cat#: 24040032). T-cells were grown and transfected in a culture medium supplemented with 10% FBS (Omega Scientific, Cat#: FB-02). When transfecting cells using mRNA-LNPs, it is typically necessary for the cell culture medium to be supplemented with 10% FBS at the time of transfection. Without this supplement, transfection efficiency will drop significantly. For mRNA-LNP transfection of cells which cannot use FBS in the culture medium, please contact us at (510) 860-4615. Alternative transfection methods are available.

## MAGE3/MHC mRNA-LNP

Ready-to-use lipid nanoparticles

### MAGE3/MHC Protein Expression in HEK293s Cells 72 Hours Post Transfection with PM-LNP-0015

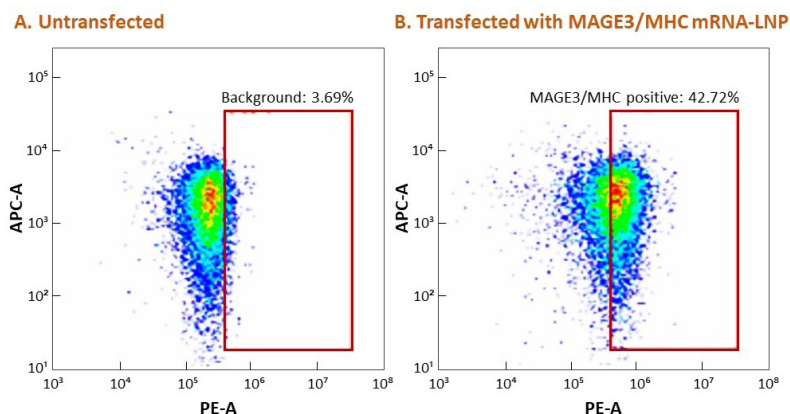


Figure 1. Flow Cytometry. PM-LNP-0015 nanoparticle-treated HEK293S cells express MAGE3/MHC protein, detected with a labeled anti-human HLA-A2 antibody (BioLegend).

### Antibody Titer

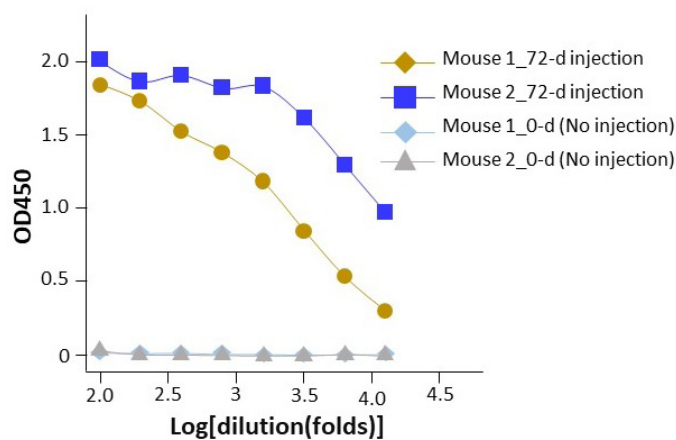


Figure 2. ELISA. Serum from mice immunized with PM-LNP-0015 nanoparticles binds to plates coated with HIA-MAGE protein, indicated by staining with a labeled anti-mouse IgG antibody.

#### Products and Services

- Monoclonal Antibody
- Bispecific Antibody
- mRNA-LNP
- Hybridoma Sequencing
- CAR-T/NK
- Stable Cell Line