

Protocol

LENTIVIRUS PRECIPITATION

OVERVIEW

The Lentivirus Precipitation Solution (cat. # VC100, ProMab Biotechnologies) is a mixture of polymer optimized for the precipitation of lentiviral particles. It provides a simple, fast and highly efficient method for concentrating lentiviral particles. In the simple protocol, you just mix your lentiviral supernatant with the Lentivirus Precipitation Solution, incubate for a short period, and spin the mixture in a standard centrifuge. You'll increase your lentivirus titer by up to 100-fold as quick as in 4 hours and obtain excellent recoveries without ultracentrifugation. The Lentivirus Precipitation Solution is a 5x solution.

PRECIPITATION PROCEDURE

- 1. Transfer the media containing lentiviral particles from plates to a sterile vessel and centrifuge the medium at 300 x g for 10 min to remove cell debris.
- 2. Filter the supernatant through 0.45 μ m filter.
- 3. Transfer filtered supernatant to a sterile vessel and add 1 volume of cold Lentivirus Precipitation Solution (4°C) to every 4 volumes of lentivirus-- containing supernatant. (Example: 5ml Lentivirus Precipitation Solution with 20ml viral supernatant).
- 4. Mix well and refrigerate 4 hours to overnight. Lentivirus-containing supernatant mixed with Lentivirus Precipitation Solution are stable for up to 4 days at 4°C.
- 5. Centrifuge mixture at 1500 × g for 30 minutes at 4°C. After centrifugation, the lentiviral particles may appear as a beige or white pellet at the bottom of the vessel.
- 6. Discard supernatant. Spin down residual solution by centrifugation at 1500 × g for 5 minutes. Remove all traces of fluid by aspiration, taking great care not to disturb the precipitated lentiviral particles in pellet.
- 7. Return the dish to cell culture incubator at 37° C with 5% CO₂.
- 8. Aliquot in cryogenic vials and store at 80°C until ready for use.

Additional Products and Services:

- Mouse Monoclonal Antibody
- Rat Monoclonal Antibody
- Rabbit Monoclonal Antibody
- Human Monoclonal Antibody
 - Polyclonal Antibody
- Antibody Sequencing
- Hybridoma Sequencing
 - CAR T-cells
 - Lentivirus production
- Cancer Stem Cells
- Specialty Cell Culture Media
- T-cell Expansion beads

Ask about our full line of CRO services to provide supplemental assistance or the entire support necessary to complete your project on time and with the data you need to move forward.

